



Inverted Research Microscope ECLIPSE Ti2

Nikon <sup>(th)</sup>  
100  
anniversary

# ECLIPSE Ti2

Inverted Research Microscope



# See More Than Be

## Leading platform for advanced imaging

The ECLIPSE Ti2 delivers an unparalleled 25mm field of view (FOV) that revolutionizes the way you see. With this incredible FOV, the Ti2 maximizes the sensor area of large-format CMOS cameras without making compromises, and significantly improves data throughput.

The Ti2's exceptionally stable, drift-free platform is designed to meet the demands of super-resolution imaging while its unique hardware-triggering capabilities enhance even the most challenging, high-speed imaging applications. Furthermore, the Ti2's unique, intelligent functions guide users through imaging workflows by gathering data from internal sensors, eliminating the possibility of user errors. In addition, the status of each sensor is automatically recorded during acquisition, providing quality control for imaging experiments and enhancing data reproducibility.

In combination with Nikon's powerful acquisition and analysis software, NIS-Elements, the Ti2 is a total innovation in imaging.

## Contents

---

Groundbreaking FOV	P4	Intuitive operation	P12
Unsurpassed Nikon optics	P6	System expandability	P14
Focus perfected	P8	NIS-Elements	P16
Assist Guide	P10	System diagrams	P18

fore



**ECLIPSE  
Ti2-E**





**ECLIPSE  
Ti2-A**



Motorized and intelligent model for advanced imaging applications. Compatible with PFS, auto correction collar, and external phase contrast system. The base of choice for live-cell imaging, high-content applications, confocal and super-resolution.

Manual model with imaging capability for laser applications. Intelligent features provide interactive guidance through imaging workflows and automatically detect microscope status.

# Groundbreaking FOV

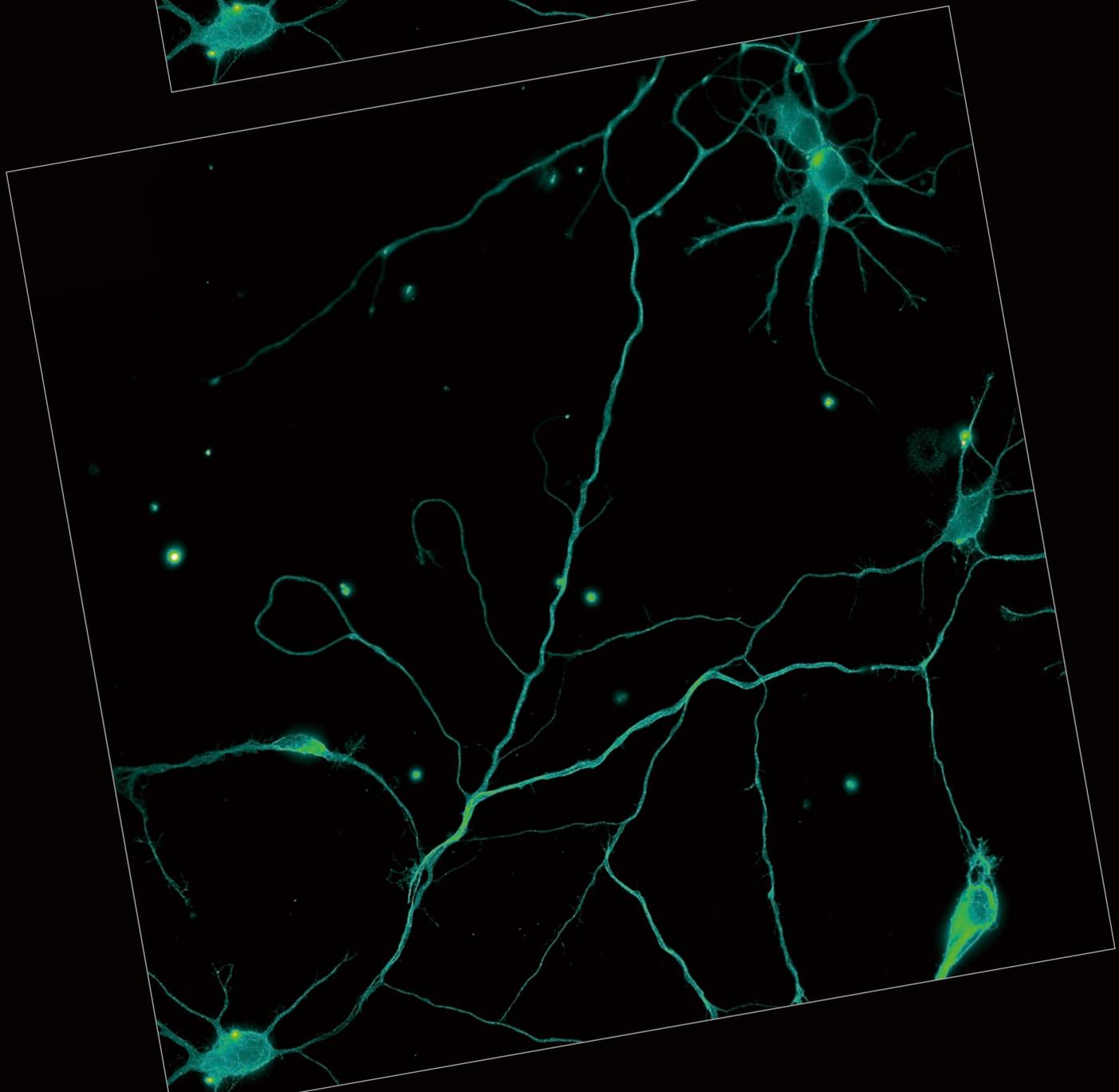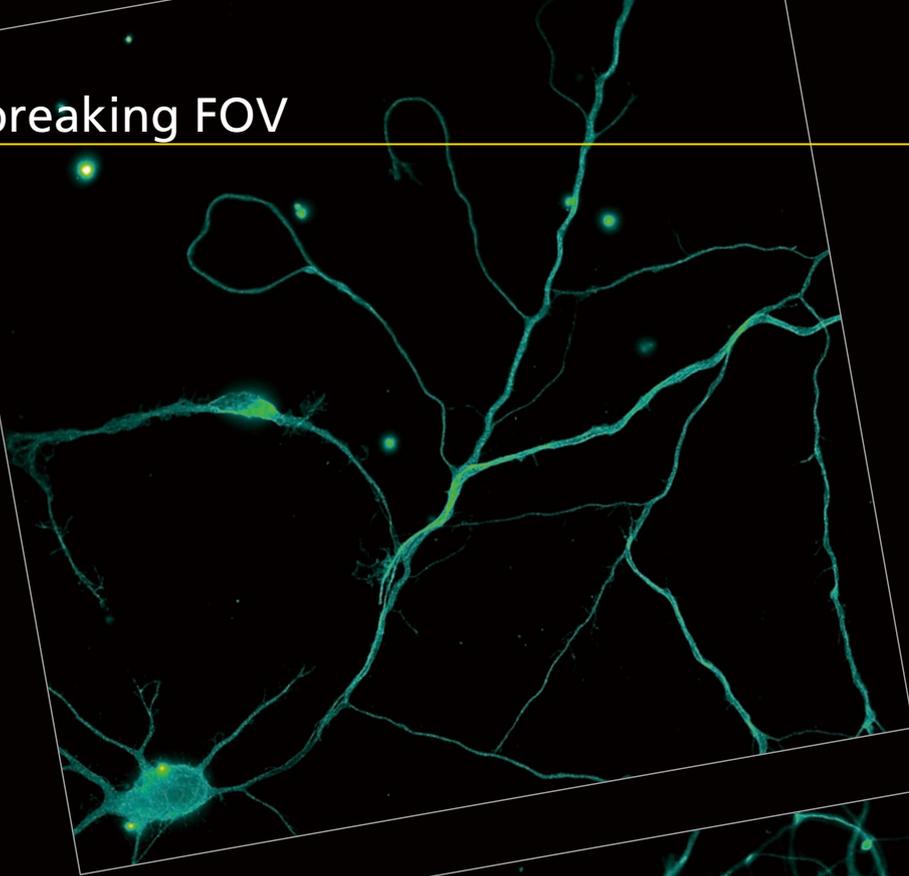


As research trends evolve towards large-scale, systems-level approaches, there is an increasing demand for faster data acquisition and higher throughput capabilities. Development of large-format camera sensors and improvements in the data processing capabilities of PCs have facilitated such research trends. The Ti2, with its unprecedented 25mm field of view, provides the next level of scalability, enabling researchers to truly maximize the utility of large-format detectors and future-proof their core imaging platform as camera technologies continue to develop at a rapid pace.

## Bright illumination over a wide area

High-power LEDs deliver bright illumination across the Ti2's large field of view, ensuring clear, consistent results from demanding applications such as high-magnification DIC. Incorporation of a fly-eye lens design provides uniform illumination from edge to edge for quantitative high-speed imaging and seamless tiling of images in stitching applications.



High-power LED illuminator

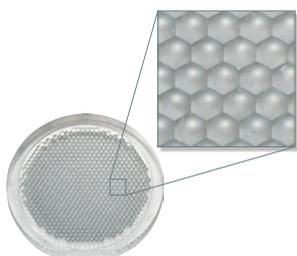

Built-in fly-eye lens

A compact epi-fluorescence illuminator designed for large FOV imaging is equipped with a quartz fly-eye lens and provides high transmittance across a broad spectrum, including UV. Large diameter fluorescence filters with hard coatings deliver large FOV images with a high signal-to-noise ratio.



Large FOV epi-fluorescence illuminator



Large diameter fluorescence filter cubes

## Large diameter observation optics

The diameter of the observation light path has been enlarged in order to achieve a field number of 25 at the imaging port. The resulting large FOV is capable of capturing approximately double the area of conventional optics, enabling users to gain maximum performance from large-format sensors such as CMOS detectors.



Enlarged tube lens



Imaging port with large 25 field number

## Objectives for large FOV imaging

Objectives with superior image flatness ensure high quality images from edge to edge. Utilizing the maximum potential of the OFN25 objective significantly accelerates data collection.



## Cameras for large-volume data acquisition

The DS-Qi2 high-sensitivity monochrome camera and DS-Ri2 high-speed color camera are equipped with large 36.0 x 23.9 mm, 16.25 megapixel CMOS image sensors, enabling maximum performance with the Ti2's large 25mm FOV.



DS-SLR camera technology optimized for microscopy



DS-Qi2








DS-Ri2

Image on page 4

Neuron culture stained for microtubules (Alexa Fluor 488), captured with CFI Plan Apo lambda 60x objective and DS-Qi2 camera. Conventional FOV on top and new Ti2 FOV on bottom.

Photo courtesy of Josh Rappoport, Nikon Imaging Center, Northwestern Univ.; Sample courtesy of S. Kemal, B. Wang, and R. Vassar, Northwestern Univ.

# Unsurpassed Nikon optics

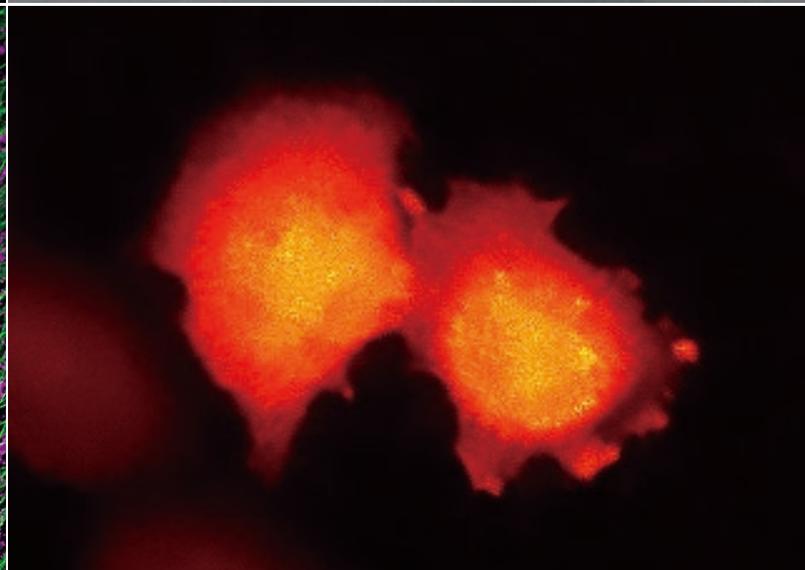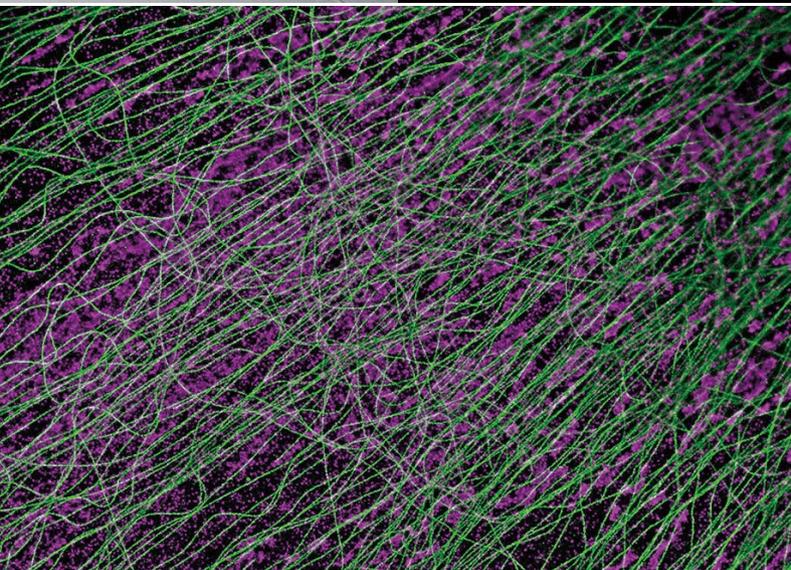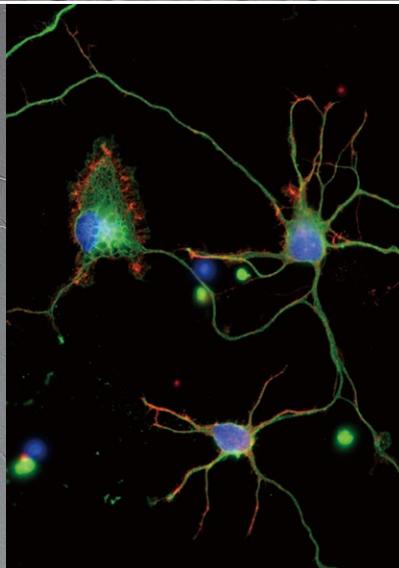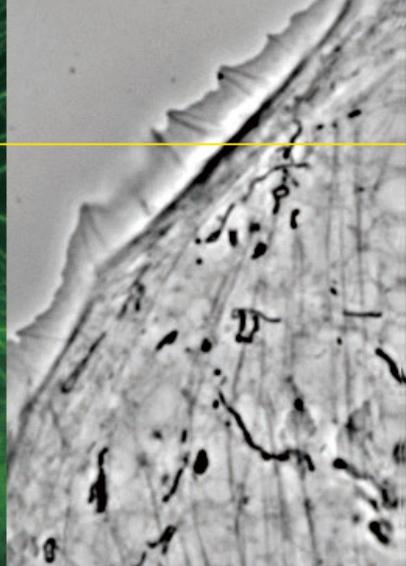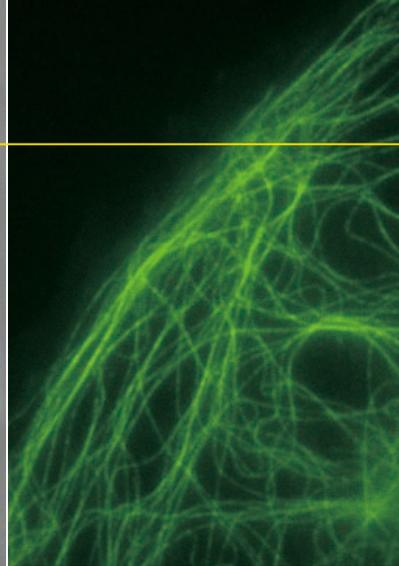

Apodized phase contrast image:  
BSC-1 cells captured with CFI S Plan Fluor ELWD ADM 40x objective

DIC and epi-fluorescence images:  
25mm FOV image of neurons (DAPI, Alexa Fluor 488, Rhodamine-Phalloidin), captured with CFI Plan Apo lambda 60x objective and DS-Qi2 camera  
Photo courtesy of Josh Rappoport, Nikon Imaging Center, Northwestern Univ.;  
Sample courtesy of S. Kemal, B. Wang, and R. Vassar, Northwestern Univ.

Super-resolution image (DNA PAINT):  
CV-1 cells expressing  $\alpha$ -tubulin (green) and TOMM-20 (magenta) captured with CFI Apo TIRF 100x Oil objective.

Epi-fluorescence and external phase contrast images:  
PTK-1 cells labeled with GFP-alpha-tubulin captured with CFI Apo TIRF 100x Oil objective  
Photo courtesy of Alexey Khodjakov, Ph.D Research Scientist VI / Professor, Wadsworth Center


NAMC image:  
Mouse embryos, captured with CFI S Plan Fluor ELWD NAMC 20x objective

Luminescence image:  
HeLa cells expressing BRET-based calcium indicator protein, Nano-lantern ( $Ca^{2+}$ ).  
Sample courtesy of Prof. Takeharu Nagai, The Institute of Scientific and Industrial Research, Osaka University

Nikon's high-precision CFI60 infinity optics, designed for use with a variety of sophisticated observation methods, are highly regarded by researchers for their superb optical performance and solid reliability.

### Apodized phase contrast

Nikon's unique apodized phase contrast objectives with selective amplitude filters dramatically increase contrast and reduce halo artifacts to provide detailed high-definition images.

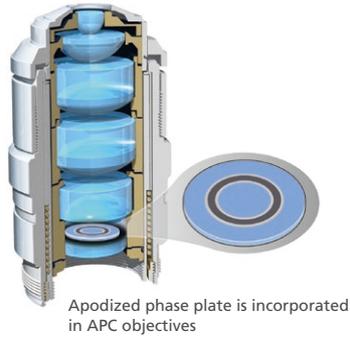

### DIC (Differential Interference Contrast)

Nikon's highly-regarded DIC optics provide uniformly clear and detailed images with high resolution and contrast throughout the magnification range. DIC prisms are individually tailored for each objective lens to provide the highest-quality DIC images for every sample.



DIC prisms matched to individual objectives are mounted in the nosepiece

### Auto Correction Collar (Ti2-E)


Changes in sample thickness, cover glass thickness, refractive index distribution in the sample, and temperature can lead to spherical aberration and image deterioration. The highest quality objectives are often equipped with correction collars to compensate for these changes, and precise positioning of the collar is critical in achieving high resolution, high contrast images. This new auto correction collar utilizes a harmonic drive and automatic correction algorithm that enable users to easily achieve precise collar adjustment to achieve optimum objective performance every time.



Harmonic drive mechanism for high-precision control of correction collar movement


### External phase contrast (Ti2-E)

The motorized external phase-contrast system enables users to combine phase contrast with epi-fluorescence imaging without compromising fluorescent light transmission by bypassing the need to use phase-contrast objectives. For example, very high NA, liquid immersion objectives can be used for phase-contrast imaging. Using this external phase contrast system, users can easily combine phase contrast with other imaging modalities, including weak-fluorescence imaging such as TIRF and laser tweezer applications.

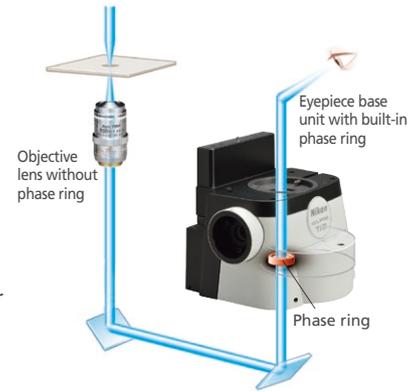


### NAMC (Nikon Advanced Modulation Contrast)

This is a plastic-compatible, high-contrast imaging technique for unstained, transparent samples such as oocytes. NAMC provides pseudo-three-dimensional images with a shadow-cast appearance. The direction of contrast can be easily adjusted for each sample.




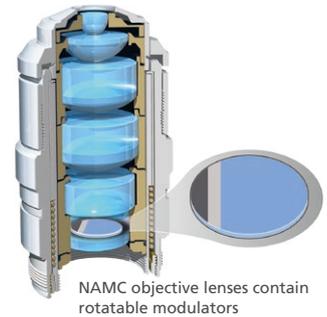


### Epi-fluorescence

The  $\lambda$  series objectives, utilizing Nikon's proprietary Nano Crystal Coat technology, are perfect for demanding, low-signal, multi-channel fluorescence imaging that requires high transmission and aberration correction over a wide wavelength range. Combined with new fluorescence filter cubes that offer improved fluorescence detection and stray light countermeasures such as the Noise Terminator, the  $\lambda$  series objectives demonstrate their power in weak signal observations such as single-molecule imaging and even luminescence-based applications.

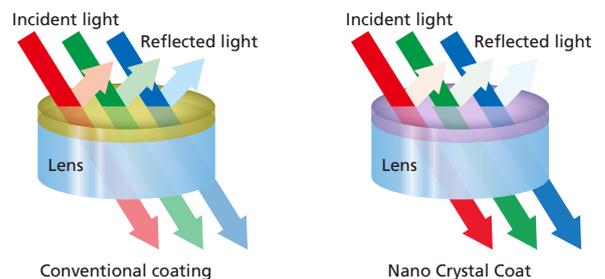


Focus perfected

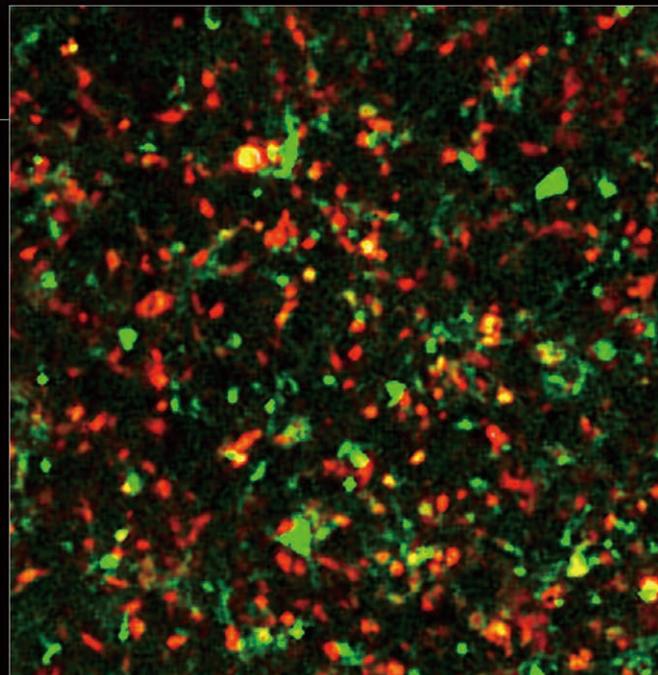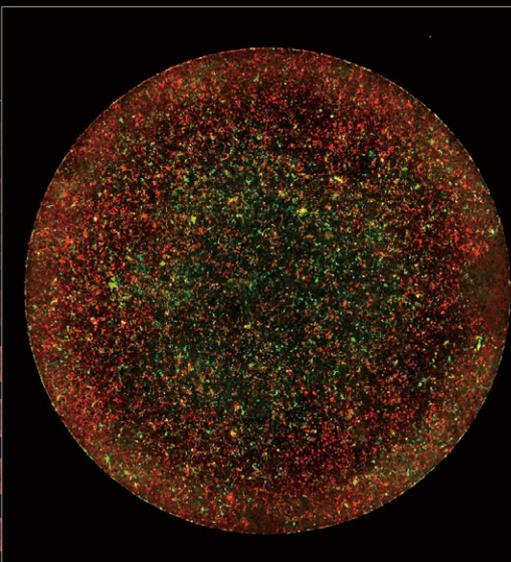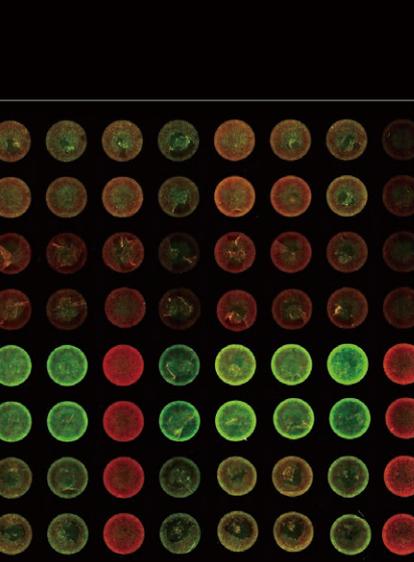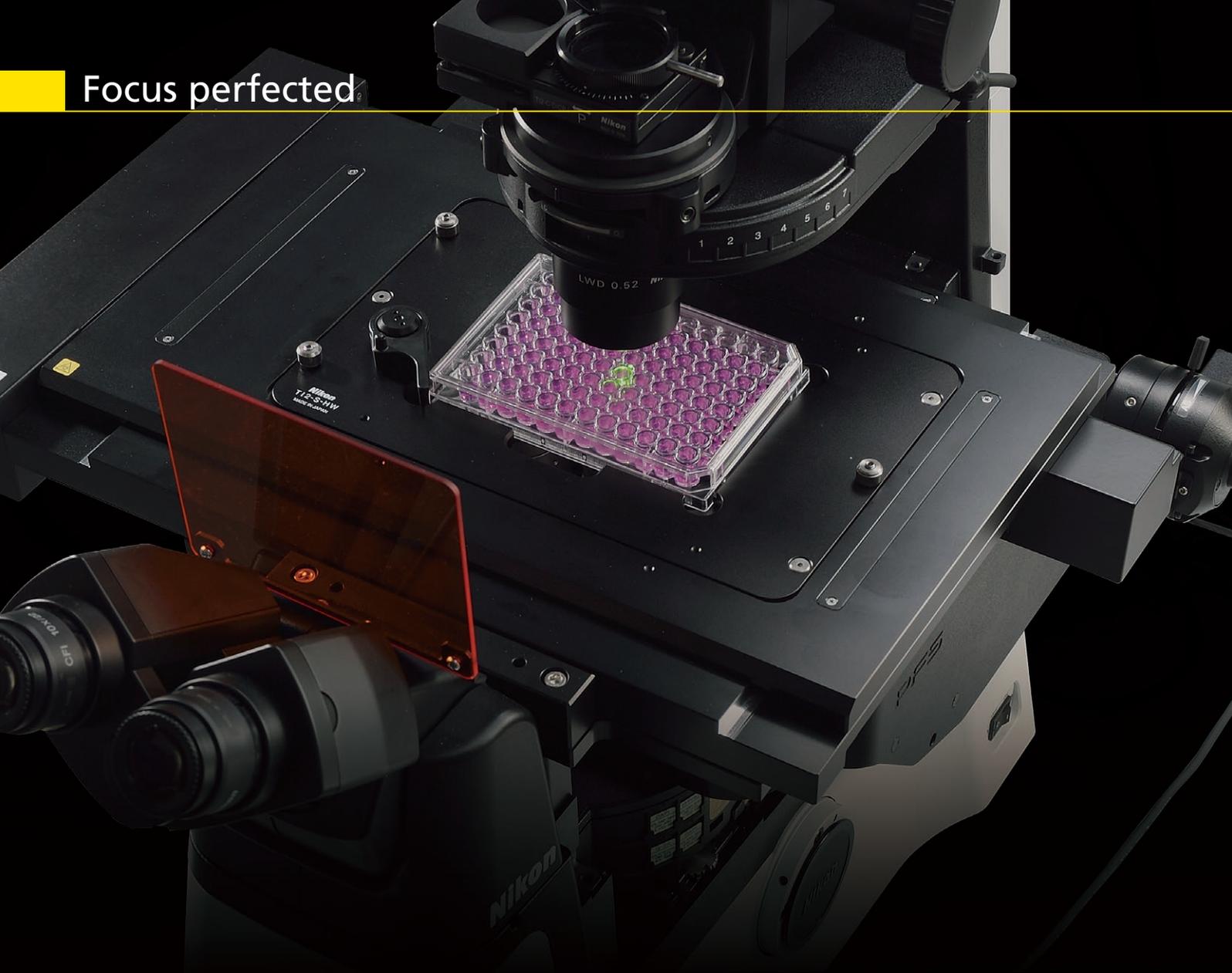

Even the slightest change in temperature and vibrations in the imaging environment can greatly impact focus stability. The Ti2 eliminates focus drift using both static and dynamic measures to enable faithful visualization of the nanoscopic and microscopic world during long time-lapse experiments.

## Mechanically redesigned for ultra-high stability (Ti2-E)

In order to improve the focusing stability, both Z-drive and PFS autofocusing mechanisms have been completely re-designed. The new Z-focusing mechanism is smaller and positioned adjacent to the nosepiece to minimize vibrations. It remains adjacent to the nosepiece even in an expanded (staged-up) configuration, ensuring stability for all applications.

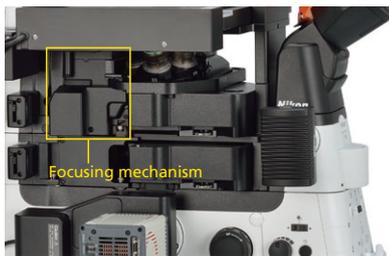

High stability Z-focusing mechanism remains adjacent to the nosepiece even in expanded configurations

The detector portion of the Perfect Focus System (PFS) has been detached from the nosepiece in order to reduce mechanical load on the objective nosepiece. This new design also minimizes heat transfer, which contributes to a more stable imaging environment. Towards this end, the power consumption of the Z-drive motor has also been reduced.


Combined, these mechanical redesigns result in an ultra-stable imaging platform, perfectly suited for single-molecule imaging and super-resolution applications.




## New generation of autofocusing with PFS: Simply perfect (Ti2-E)

The newest generation of the Perfect Focus System (PFS) automatically corrects focus drift caused by temperature changes and mechanical vibrations that can be caused by a variety of factors, including the addition of reagents to the sample and multi-position imaging.

PFS detects and tracks the position of a reference plane (e.g. coverslip surface in the case of immersion objectives) in real time to maintain focus. Unique optical-offset technology allows users to maintain focus at an arbitrary position offset from the reference plane. The user can simply focus on the desired plane and then engage PFS. PFS then automatically and continuously maintains focus by means of a built-in linear encoder and high-speed feedback mechanism, providing highly-reliable images even during long-term, complex imaging tasks.

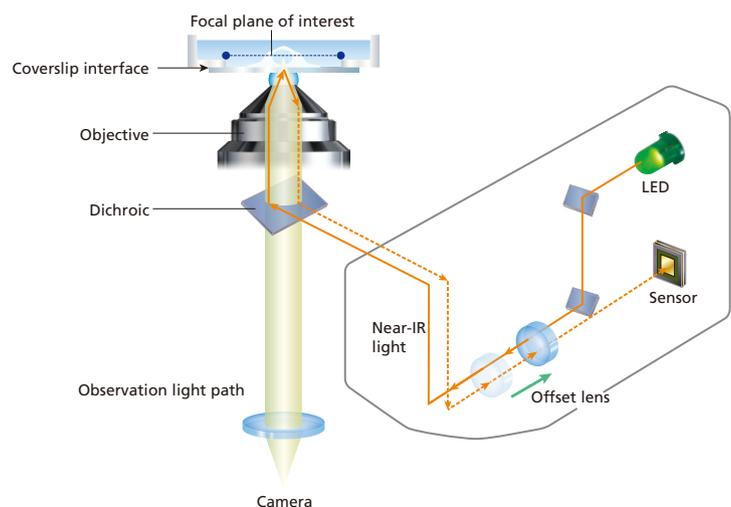

PFS is compatible with a wide range of applications, from routine experiments involving plastic culture dishes to single-molecule imaging and multi-photon imaging. It is also compatible with a wide range of wavelengths, from ultraviolet to infrared, meaning it can be used for multi-photon and optical tweezer applications.

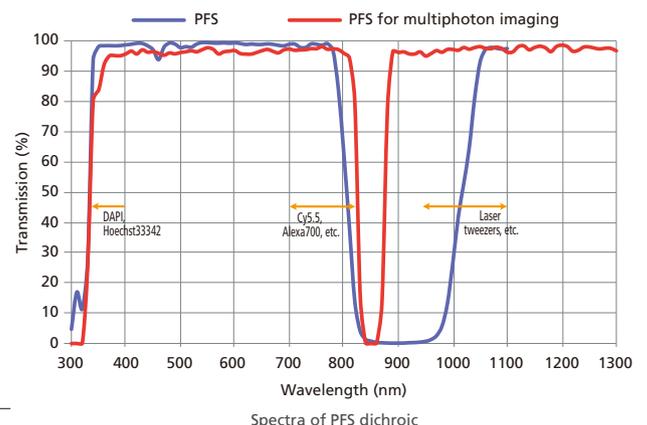

Images on page 8  
 Stitched large image of an entire 96-well plate using PFS and a CFI Plan Apo lambda 4x objective; each well contains neuronal cultures expressing red and green fluorescent proteins.  
 Sample courtesy of Jeanette Osterloh and Steve Finkbeiner, Gladstone Institutes, UCSF



It is no longer necessary to memorize complex microscope alignment and operation procedures. The Ti2 integrates data from sensors to guide you through these steps, eliminating user error and enabling researchers to concentrate on their data.

### Continuous display of microscope status (Ti2-E/A)

A collection of built-in sensors detects and relays status information for a variety of components in the microscope. All of the status information is recorded in the metadata when you acquire images with a computer, so you can easily recall acquisition conditions and/or check for configuration errors.

In addition, a built-in internal camera allows users to view the back aperture, facilitating confirmation of phase ring alignment and extinction cross in DIC. It also provides a laser-safe method for aligning lasers for applications such as TIRF.

Microscope status can be viewed on a tablet and also determined based on status lights on the front of the microscope, enabling status determination in a dark room.




Status lights



Built-in sensors detect the status of microscope components


### Guidance for operational procedures (Ti2-E/A)

The Ti2's Assist Guide function provides interactive step-by-step guidance for microscope operation. The Assist Guide can be viewed on a tablet or PC, and integrates real time data from built-in sensors and an internal camera. The Assist Guide is designed to help users through alignment procedures for both experiment setup and troubleshooting.

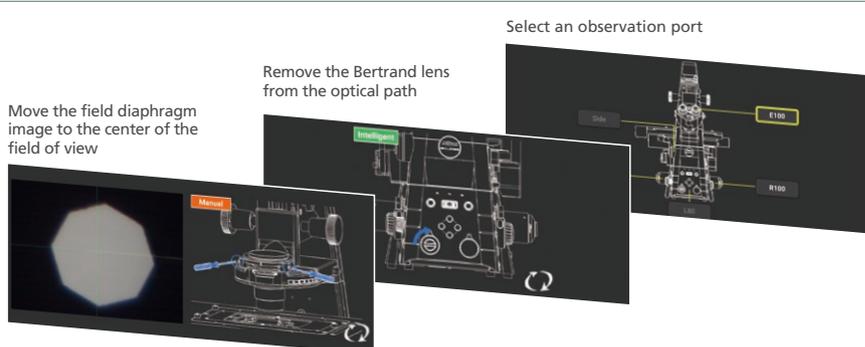

### Automatically detect errors (Ti2-E/A)

The Check Mode allows users to easily confirm, on either a tablet or PC that all the correct microscope components are in place for their chosen observation method. This capability eliminates time and effort normally required for troubleshooting when the desired observation method is not achieved. This functionality is particularly advantageous when multiple users are involved, each with the potential to make unexpected changes to the microscope settings. Custom check procedures can also be pre-programmed.

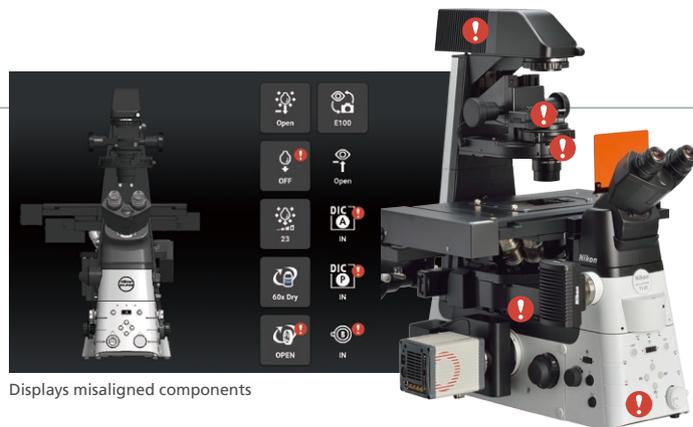


Displays misaligned components



The Ti2 has been completely redesigned, from the overall body design to the selection and placement of every button and switch, for the ultimate in user experience. The controls are easy to use even in the dark, where the majority of imaging experiments are conducted. The Ti2 provides an intuitive and effortless user interface so researchers can focus on the data and not on microscope control.

### Thoughtfully designed layout for microscope control (Ti2-E/A)

The placement of all of the buttons and switches are based on the type of illumination they control. Buttons that control diascopic observation are positioned on the left side of the microscope and those that control epi-fluorescence observation are on the right side. Buttons that control common operations are on the front panel. This use of zoning provides an easy-to-remember layout, a desirable feature when operating the microscope in a dark room.

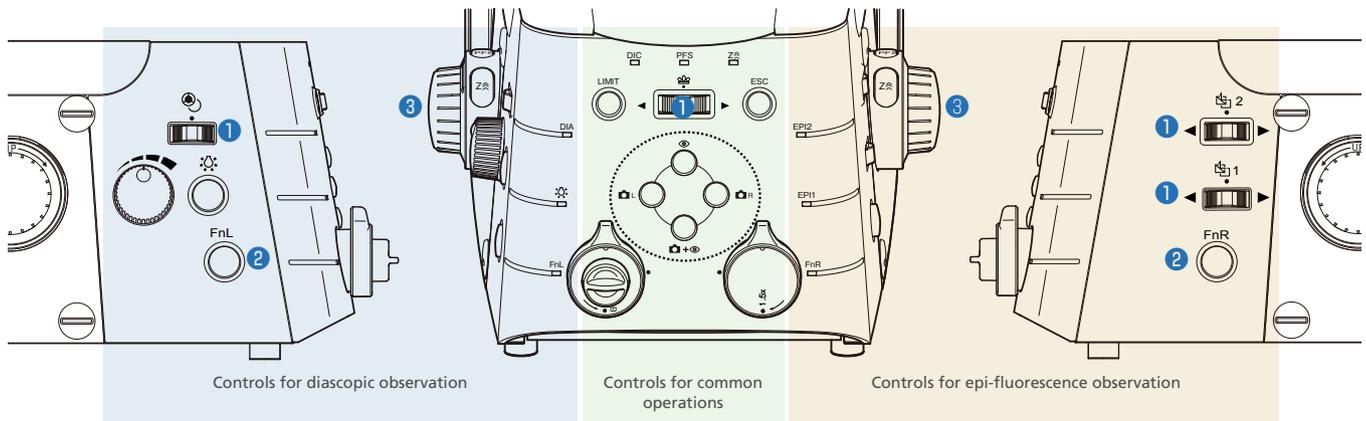

#### 1 Shuttle switch (Ti2-E)

Shuttle switches have been incorporated into the design to control devices such as the fluorescence filter turret and objective nosepiece. These types of switches emulate the feel of manually rotating these devices, for intuitive control. Additional functionality can be incorporated into these shuttle switches so that a single switch can operate multiple related devices. For example, the shuttle switch for the fluorescence filter turret not only rotates the turret but also opens and closes the fluorescence shutter when the user presses the switch. It is also possible to program these switches to operate a barrier filter wheel and the external phase contrast unit.

#### 2 Programmable Function button (Ti2-E/A)

Conveniently located Function buttons allow customization of the user interface. Users can select from more than 100 functions, including control of motorized devices such as shutters and even signal output to external devices via the I/O port for triggered acquisition. Mode functions, which enable instant changing of observation methods by storing the settings of each motorized device, can also be assigned to these buttons.

#### 3 Focusing knob (Ti2-E)

A focus acceleration button and a PFS engagement button are provided adjacent to the focusing knobs. The two buttons are easily identified by touch because of their different shapes. Focusing speed is automatically adjusted for the objective in use, enabling stress-free operation by maintaining an ideal focusing speed.

### Intuitive control with joystick and tablet (Ti2-E)

The Ti2 joystick not only controls stage movement, but the majority of motorized functions on the microscope, including PFS activity. It can display XYZ coordinates and the status of microscope components, providing an effective means for the user to remotely control the microscope. Motorized functions of the Ti2 can also be controlled from a tablet, connected by wireless LAN to the microscope, providing a versatile graphical interface for microscope control.



## System expandability



### Stage up kit

Allows expansion of the infinity space for incorporation of additional devices such as a second fluorescence filter turret, barrier filter wheel, back port unit, and LAPP modules.



### Epi-fluorescence filter turret

Compatible with large FOV. A manual intelligent model and high-speed motorized models with motorized shutter are available.



### Motorized barrier filter wheel

Provides high-speed filter switching (50msec. switch time between adjacent positions) and is compatible with large FOV. It can also be mounted in the infinity space, under the fluorescence filter turret, when the stage is raised.



### Motorized stage

Control methods are optimized for fast stage operation. A stage with a built-in encoder and a piezo stage for Z-stacking are available. Magnetic specimen holders ensure safe operation.



### Nosepiece

Manual, intelligent and motorized nosepieces, PFS-specific nosepieces including models for auto correction collar and multiphoton imaging are available. The design of these nosepieces prevents internal leakage of immersion liquids and sample media.



### Controller for Ti2-E

Allows the Ti2-E to communicate with external devices and enables high-speed triggered acquisitions. I/O ports can be added by connecting an extension box.



### LAPP system

Various illumination modules, such as TIRF and photostimulation modules, can be flexibly combined to customize your imaging system. Up to five modules can be mounted simultaneously.



### Field stop sliders

Two different rectangular-shaped apertures and one round aperture model are available. The rectangular models prevent excitation outside of the imaging area and unintentional photobleaching of samples. They can be removed for ultra-wide FOV imaging.



### Condenser turret

Three types are available: manual, intelligent, and a motorized type that automatically switches between seven positions to accommodate different imaging modes.



### Stage top incubator

The STX series precisely maintains temperature at 37.0°C and humidity at more than 95%, and controls CO<sub>2</sub> to enable culturing of cells for more than 1 week.

Manufactured by Tokai Hit Co., Ltd.



### Manual stage

A long travel stroke allows observation of the entire area of a well plate. The travel area can be restricted to ensure safe operation. Various specimen holders are available.



### Confocal microscopes

The A1+ high-resolution/A1R+ ultra-fast high-resolution confocal system and C2+ personal confocal system. Multiphoton versions are also available for the A1 confocal series.



### Oil hydraulic micromanipulator

The MTK-1-N4 achieves simple, easy pipette alignment and smooth pipette movement by means of a four-axis hydraulic joystick.

Manufactured by Narishige Co., Ltd.



### Epi-FL LED lightsource

Stable, high-intensity light source for maintenance-free, long-life illumination. The HG fiber illuminator "Intensilight" is also available.



### Super Resolution Microscopes

N-SIM, which achieves live-cell imaging with double the resolution of conventional optical microscopes, and N-STORM, which achieves ten times the resolution of conventional optical microscopes for molecular-level observation, are available.

# NIS Elements


## A unified acquisition and analysis software platform

Nikon's unified software platform, NIS-Elements, provides acquisition control for basic to advanced imaging systems as well as powerful analysis tools and stunning display capabilities. NIS-Elements can be streamlined for simple, turnkey use and expanded for fully customized, complex experiments such as conditional workflows. NIS-Elements also offers easy-to-use, graphical programming tools such as JOBS and Illumination Sequence for customizing tasks. In addition to Nikon hardware, NIS-Elements controls devices from a vast array of manufacturers to enable the highest level of customization.

## Device Control


### Multidimensional Imaging

Optical Configuration (OC) settings memorize custom observation modes and are combined in the ND Acquisition GUI to easily create experiments combining multi-channel, multi-stage position, z-stacking, and timelapse imaging, as well as image stitching. Other functions such as photostimulation and photobleaching can also be flexibly combined.

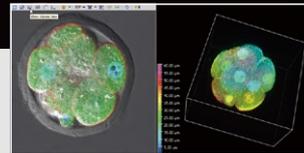


### High-speed hardware control

The Ti2's unique hardware triggering capabilities remove software callbacks during acquisition routines to maximize imaging speed. The Illumination Sequence module provides a simple, graphical interface for quickly designing and verifying complex, triggered acquisition experiments.

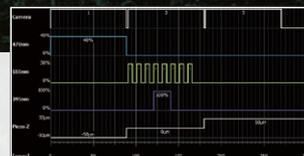


### Graphical programming for custom tasks

For complex or unique image acquisition needs, the JOBS tool provides an easy-to-use graphical interface for creating custom workflows using drag-and-drop features.

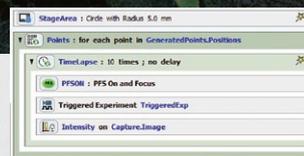

### Control of Third Party Products

NIS-Elements can natively control a vast number of devices from a variety of manufacturers including high-sensitivity cameras, piezo-devices, light sources, wheels, and National Instruments DAQ devices. This flexibility in hardware control enables custom configurations tailored to individual research needs.

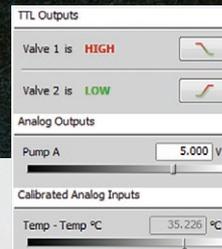


## Display & Processing

## Image Analysis

### Multi-dimensional Display

Multi-dimensional images that combine multichannel, time-lapse, Z stack, and multi-XY position imaging, as well as image stitching are intuitively displayed in a single window. Powerful volume rendering tools provide unrivalled image quality and interactivity while the Movie-Maker feature enables users to easily create stunning movies.

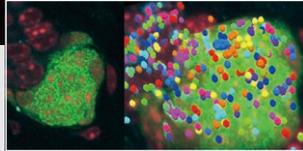


### Advanced Image Processing

Advanced filters for sharpening, smoothing, and denoising as well as real-time shading correction and image averaging for noise reduction are available. NIS-Elements also offers advanced image arithmetic tools as well as a variety of image projection options such as extended depth of focus (EDF).




### Deconvolution

Automatic and manual modes, robust algorithms for noise measurement and removal, and enhanced spherical aberration correction are provided to help actualize theoretical resolutions for even confocal images. Both 3D and 2D deconvolution options are available.

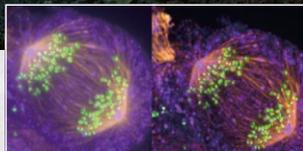


### Automatic measurement

A combination of powerful segmentation tools, morphology functions, classifiers, and an extensive list of measurement tools for 2D, 3D and timelapse datasets enable users to extract quantitative information from their data with ease. Interactive/manual measurement options are also available.

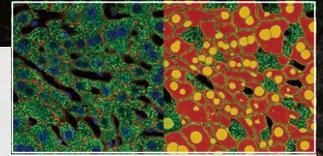

### Real-time measurement

Time measurements can be carried out in real time and visualized during acquisition. Real-time analysis results can be especially useful for optimizing experiments that rely on drug treatments or ratiometric imaging like FRET or FRAP.

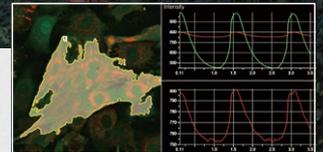


### 2D and 3D Object Tracking

NIS-Elements provides powerful tools for identifying and tracking 2D and 3D objects. Measurements include velocity, acceleration, distance, and direction. A variety of innovative display options are available for communicating and presenting tracking analysis results.



### Custom Analysis Routines




The General Analysis (GA) module provides an easy method for creating custom analysis routines that combine image processing and measurement. Routines created in GA can be saved and recalled and even combined with JOBS to create conditional acquisition workflows that rely on real-time analysis results from the integrated GA routine.

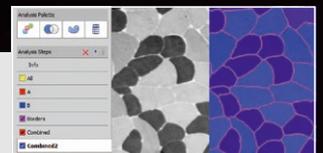


# System Diagram

Motorized accessories (with status detection function)


Accessories with status detection function

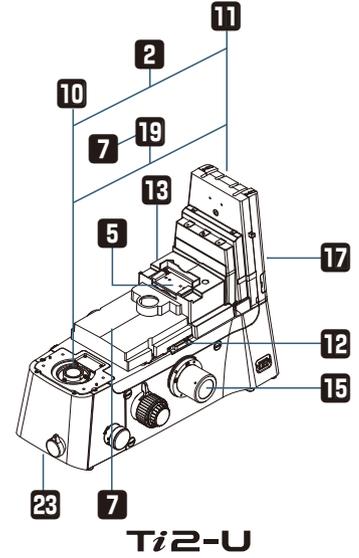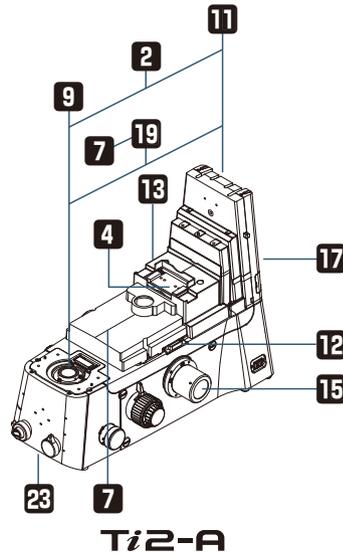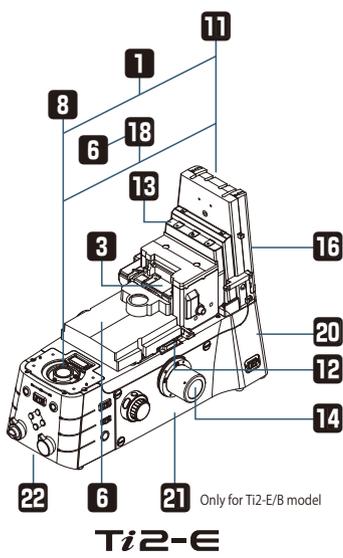


## Stages

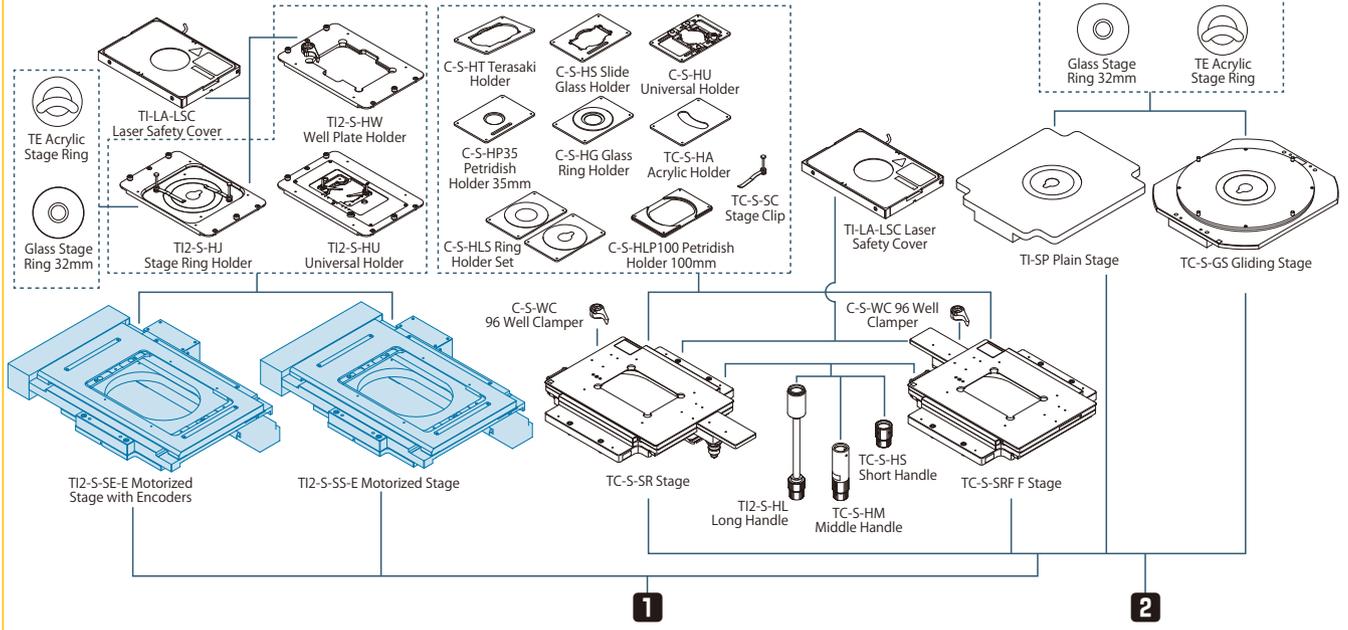


## Nosepieces

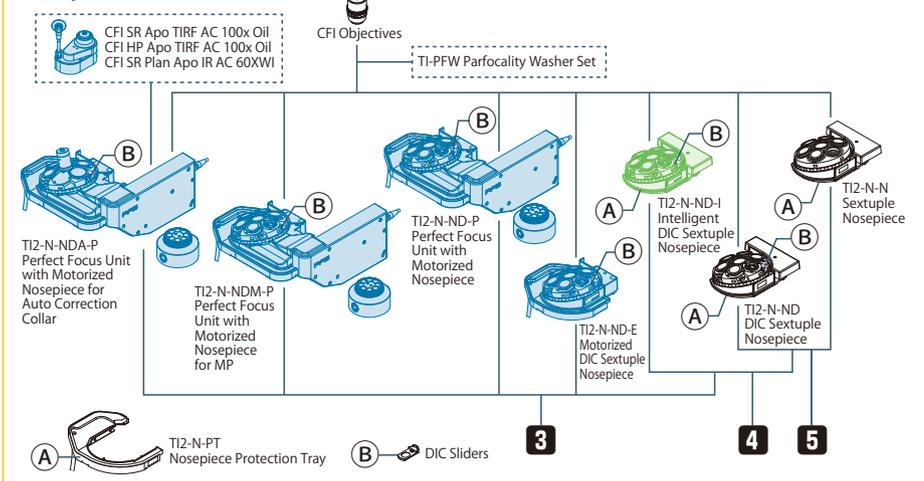


## Fluorescence Cube Turrets

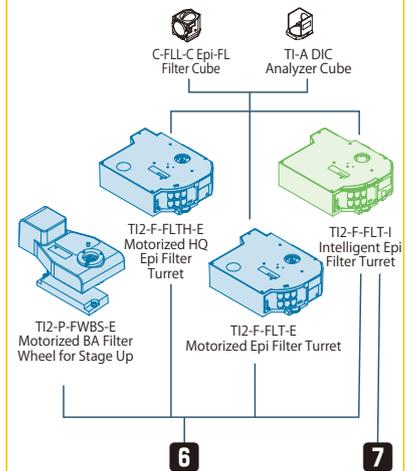


### Tubes/Eyepieces

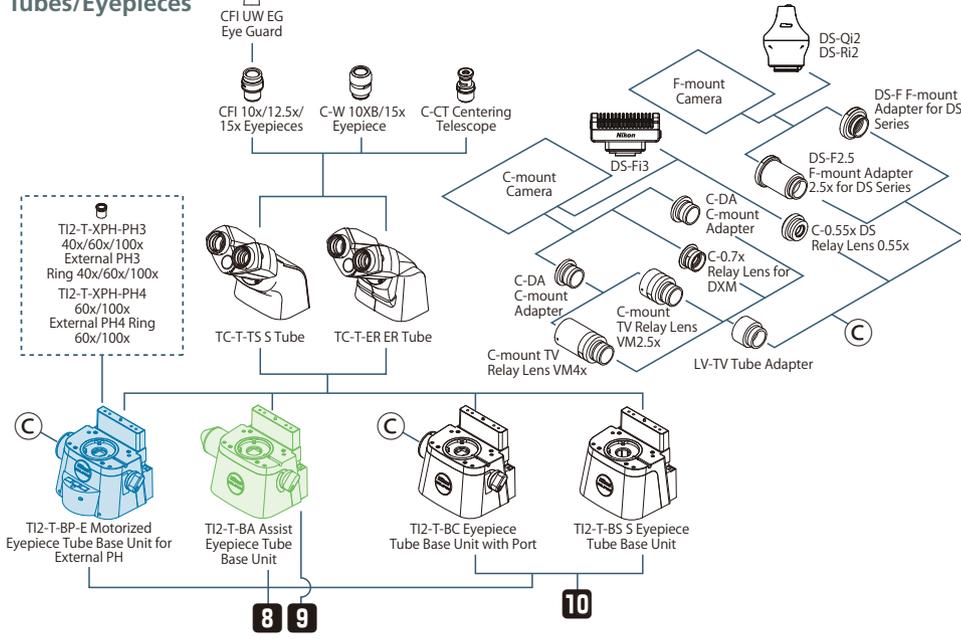


### Bottom Port

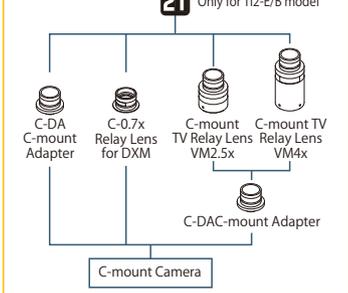


### Controllers

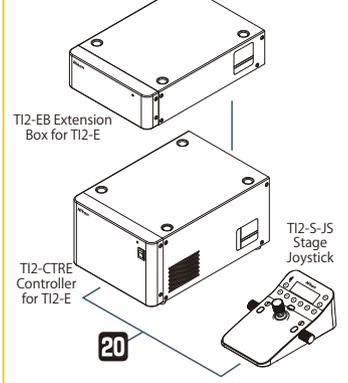


### Diascopic Illuminators

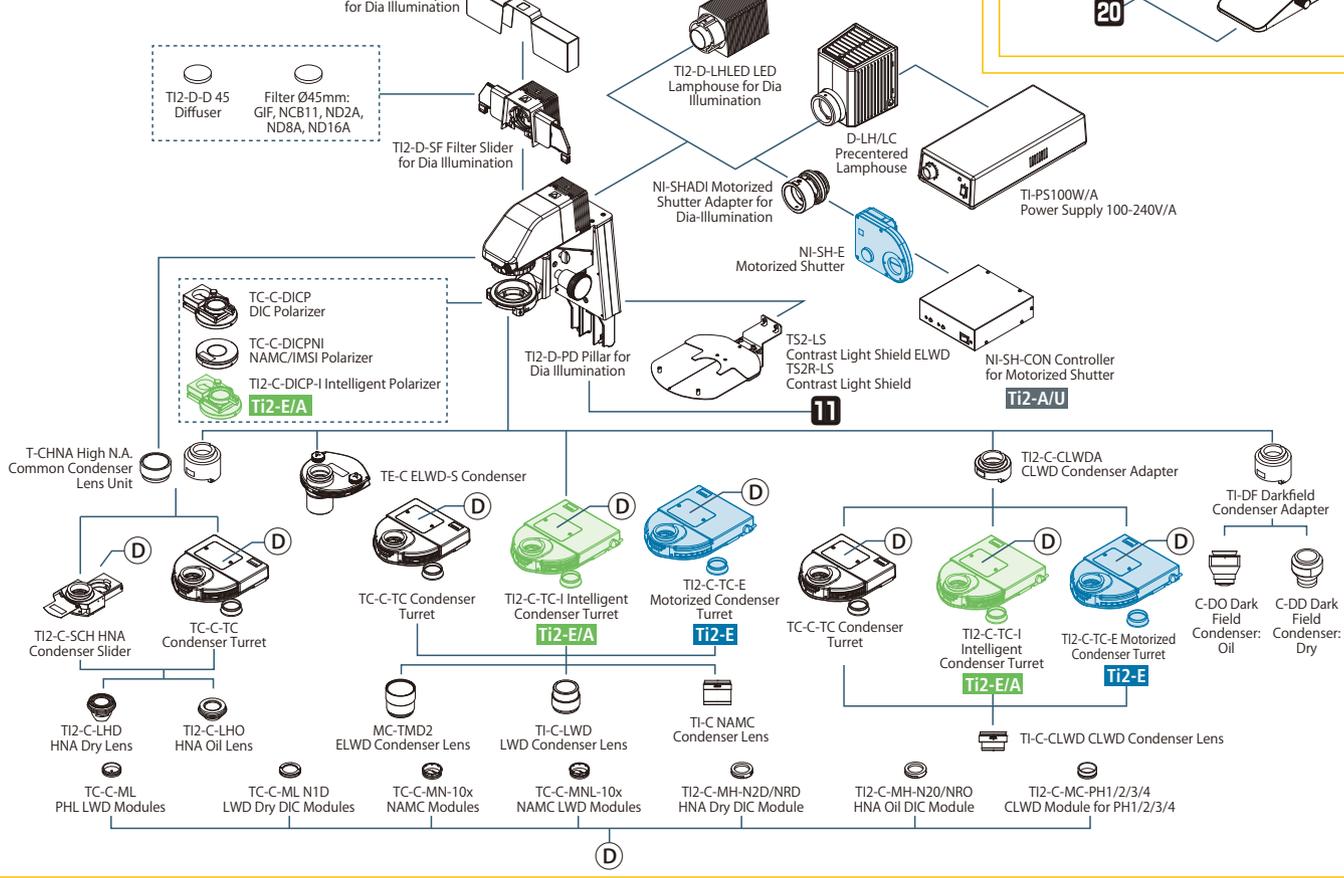



### Stage Risers

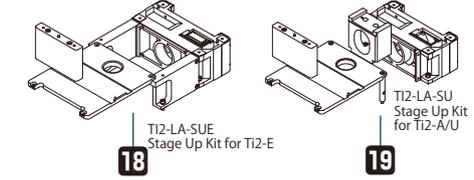

### Analyzer

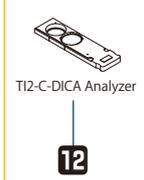

### Field Stops

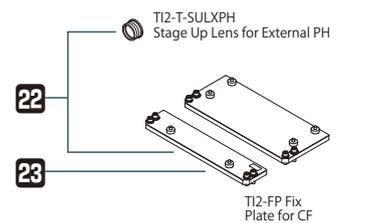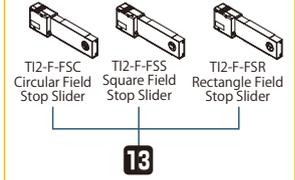


# System Diagram


Motorized accessories (with status detection function)

Accessories with status detection function

## Side Port

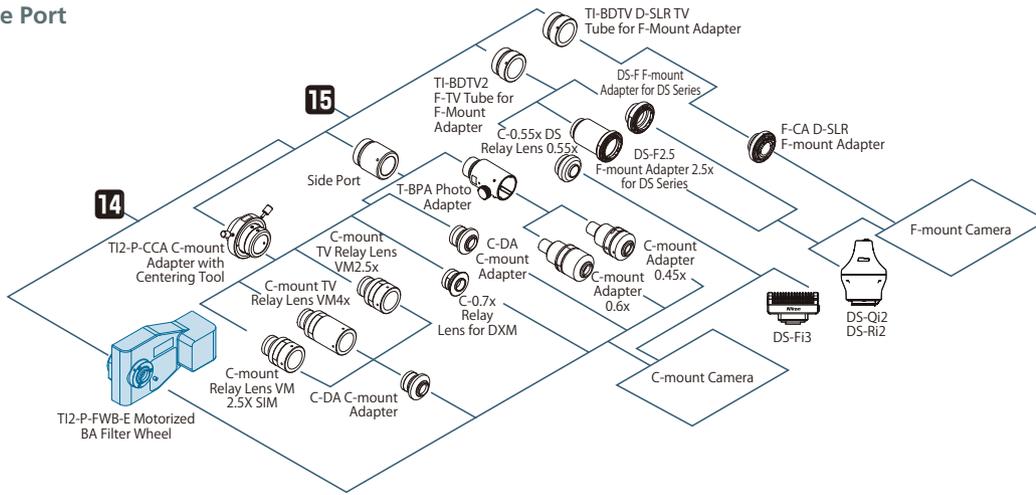

## Fluorescence Illuminators

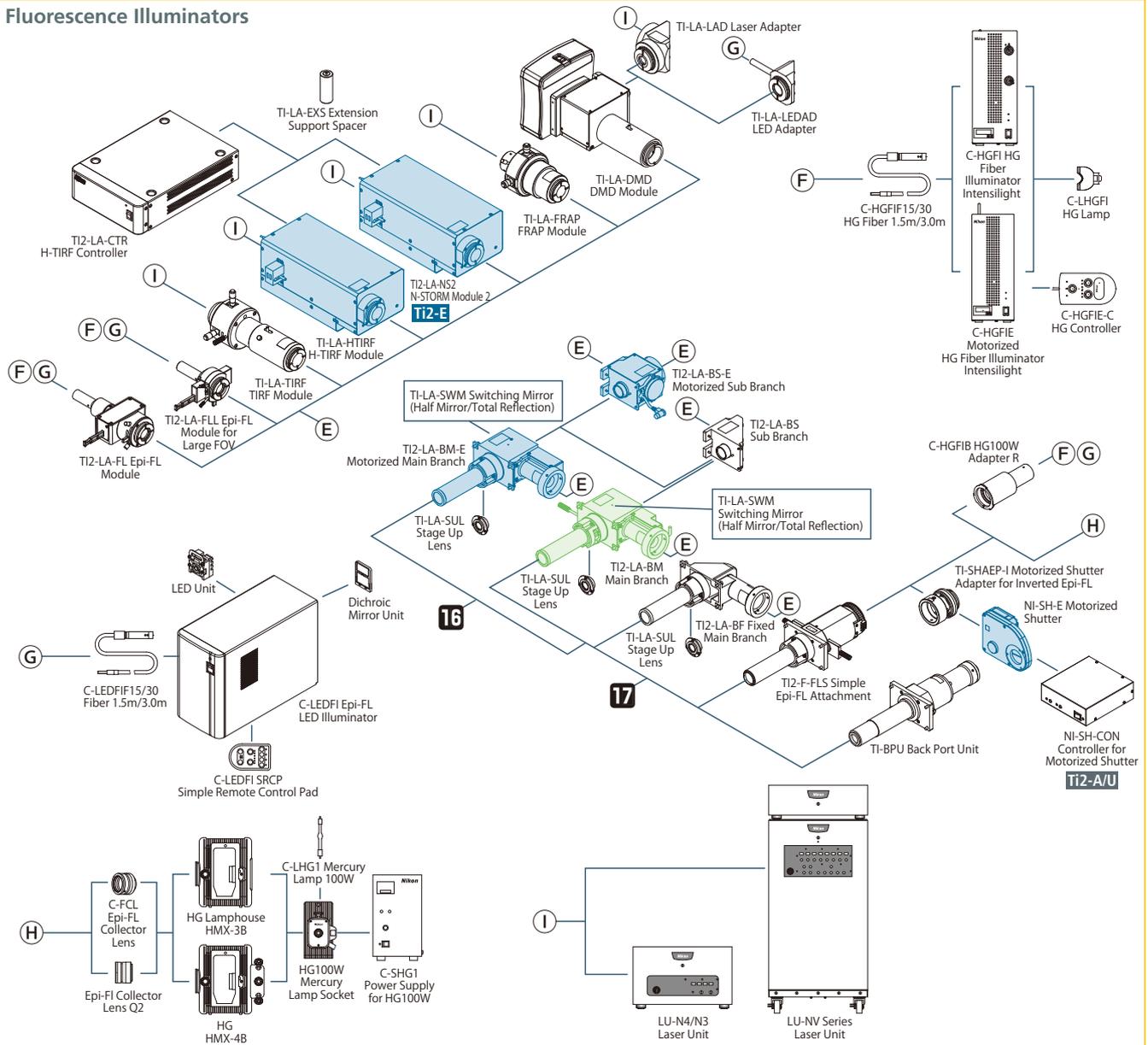

# Objectives

Type	Model	Immersion	NA	W.D. (mm)	Cover glass thickness	Correction ring	Spring loaded	DIC	Phase contrast	Fluorescence			PFS	
										UV	Visible light	NIR		
Plan Fluor	CFI Plan Fluor	4x		0.13	17.20	—					☉	☉		
		10x		0.30	16.00	0.17			○		☉	☉		✓
		20x		0.50	2.10	0.17			○		☉	☉		
		20xA MI	Oil Water Glycerin	0.75	0.51-0.35 0.51-0.34 0.49-0.33	0-0.17	✓	✓	○		☉	☉		
		40x		0.75	0.66	0.17		✓	○		☉	☉		✓
		40x Oil	Oil	1.30	0.24	0.17		✓w/stopper	○	EXT PH3-40x	☉	☉		✓
		60x		0.85	0.40-0.31	0.11-0.23	✓	✓	○		☉	☉		
		60xSH (with iris)	Oil	0.50-1.25	0.22	0.17		✓	○	EXT PH3-60x	☉	☉		
		100x Oil	Oil	1.30	0.16	0.17		✓w/stopper	○		☉	☉		✓
		100xSH (with iris)	Oil	0.50-1.30	0.16	0.17		✓	○		☉	☉		
		DL 4x		0.13	16.40	1.2				PHL	○	○		
		DLL 10x		0.30	16.00	0.17				PH1	○	○		✓
		DL 10x		0.30	15.20	1.2				PH1	○	○		✓
		DLL 20x		0.50	2.10	0.17				PH1	○	○		✓
		DLL 40x		0.75	0.66	0.17			✓	PH2	○	○		✓
		DLL 100x Oil	Oil	1.30	0.16	0.17		✓w/stopper		PH3	○	○		✓
		ADH 100x Oil	Oil	1.30	0.16	0.17		✓w/stopper		PH3	○	○		✓
S Plan Fluor	CFI S Plan Fluor	ELWD 20xC		0.45	8.2-6.9	0-2.0	✓		○		☉	☉		✓
		ELWD 40xC		0.60	3.6-2.8	0-2.0	✓		○		☉	☉		✓
		ELWD 60xC		0.70	2.6-1.8	0.1-1.3	✓		○		☉	☉		
		ELWD ADM 20xC		0.45	8.2-6.9	0-2.0	✓			PH1	○	○		✓
		ELWD ADM 40xC		0.60	3.6-2.8	0-2.0	✓			PH2	○	○		✓
		ELWD ADL 60xC		0.70	2.6-1.8	0.1-1.3	✓			PH2	○	○		
		ELWD NAMC 20xC		0.45	8.2-6.9	0-2.0	✓				○	○		
ELWD NAMC 40xC		0.60	3.6-2.8	0-2.0	✓				○	○				
Super Fluor	CFI Super Fluor	4x		0.20	15.50	—					☉ 340	☉		✓
		10x		0.50	1.20	0.17		✓	○		☉ 340	☉		✓
		20x		0.75	1.00	0.17		✓	○		☉ 340	☉		✓
		40x		0.90	0.30	0.11-0.23	✓	✓	○		☉ 340	☉		
		40x Oil	Oil	1.30	0.22	0.17		✓w/stopper	○	EXT PH3-40x	☉ 340	☉		✓
		100xSH (with iris)	Oil	0.5-1.3	0.20	0.17		✓			☉ 340	☉		
Plan Apochromat	CFI Plan Apo	λ 2x		0.10	8.50	—					●	☉	☉	
		λ 4x		0.20	20.00	—					●	☉	☉	✓
		λ 10x		0.45	4.00	0.17		✓	○		●	☉	☉	✓
		10xC Glyc	Oil Water Glycerin	0.5	5.50	0-0.17	✓				●	☉	☉	
		λ 20x		0.75	1.00	0.17		✓	○		●	☉	☉	✓
		VC 20x		0.75	1.00	0.17		✓	○		●	☉	☉	✓
		λ 40x		0.95	0.25-0.17	0.11-0.23	✓	✓	○		●	☉	☉	✓
		λ 60x		0.95	0.20-0.11	0.11-0.23	✓	✓	○		●	☉	☉	
		λ 60x Oil	Oil	1.40	0.13	0.17		✓	○	EXT PH3-60x	●	☉	☉	✓
		VC 60xA WI	Water	1.20	0.31-0.28	0.15-0.18	✓	✓	○	EXT PH3-60x	●	☉	☉	✓
		IR 60xWI	Water	1.27	0.18-0.16	0.15-0.19	✓	✓	○	EXT PH3-60x	●	☉	☉	✓
		λ 100x Oil	Oil	1.45	0.13	0.17		✓	○	EXT PH3-100x	●	☉	☉	✓
		VC 100x Oil	Oil	1.40	0.13	0.17		✓	○	EXT PH3-100x	●	☉	☉	✓
		HP VC 100x Oil	Oil	1.40	0.13	0.17		✓	○	EXT PH3-100x	●	☉	☉	✓
		λ DM 20x		0.75	1.00	0.17		✓		PH2	●	○	○	✓
		λ DM 40x		0.95	0.25-0.17	0.11-0.23	✓	✓		PH2	●	○	○	✓
		λ DM 60x		0.95	0.20-0.11	0.11-0.23	✓	✓		PH2	●	○	○	
		λ DM 60x Oil	Oil	1.40	0.13	0.17		✓		PH3	●	○	○	✓
		λ DM 100x Oil	Oil	1.45	0.13	0.17		✓		PH3	●	○	○	✓
		SR IR 60xWI*	Water	1.27	0.18-0.16	0.15-0.19	✓		○	EXT PH3-60x	●	○	☉	✓
Apochromat	CFI Apo	LWD 20xWI λS	Water	0.95	0.95	0.11-0.23	✓		○		○	○	✓	
		40xWI λS	Water	1.25	0.20-0.16	0.15-0.19	✓	✓	○	EXT PH3-40x	☉	☉	✓	
		LWD 40xWI λS	Water	1.15	0.60	0.15-0.19	✓	✓	○	EXT PH3-40x	○	☉	✓	
		60x Oil λS	Oil	1.40	0.14	0.17		✓	○	EXT PH3-60x	☉	☉	✓	
		TIRF 60x Oil	Oil	1.49	0.12	0.13-0.19 (23°C) 0.15-0.21 (37°C)	✓		○	EXT PH4-60x	●	☉		✓
		TIRF 100x Oil	Oil	1.49	0.12	0.13-0.19 (23°C) 0.14-0.20 (37°C)	✓		○	EXT PH4-100x	●	☉		✓
		HP TIRF 100x Oil*	Oil	1.49	0.12	0.13-0.19 (23°C) 0.14-0.20 (37°C)	✓		○	EXT PH4-100x	●	☉		✓
		SR TIRF 100x Oil*	Oil	1.49	0.12	0.13-0.19 (23°C) 0.14-0.20 (37°C)	✓		○	EXT PH4-100x	●	☉		✓

☉ : recommended for best results

○ : suitable

● : possible with visible light that has a longer wavelength than excitation light used for DAPI

EXT : external phase contrast modules

340 : high transmittance with an ultraviolet wavelength range of up to 340nm

\* Objectives for Auto Correction Collar are available.

# Specifications

		ECLIPSE Ti2-E/Ti2-E/B*1	ECLIPSE Ti2-A	ECLIPSE Ti2-U
Main body	Optical system	Infinity-corrected CFI60		
	Field number*2	22 with C-mount, 25 with F-mount		
	Intermediate Magnification switching	Manual switching of 1.0x/1.5x (exchangeable from 1.5x to 2.0x)		
		Status detection	—	
	Bertrand lens	Manual in/out, and manual focus, Status detection		
	Output port	4 Motorized positions Eyepiece 100%, left 100%, right 100%, eyepiece 20%/left 80% (Ti2-E/B: bottom 100%)	4 Manual positions Eyepiece 100%, left 100%, right 100%, option (to eyepiece 20%/left 80% or eyepiece 20%/right 80%)	
		Can add ports by use of back port unit and/or choice of tube base unit*3		
Focusing unit	Motorized drive, Coarse/fine focus changeover, 10mm stroke, Minimum increments: 0.01µm, 0.02µm (with encoder control)	Manual drive, Coarse/fine focusing knob, 10mm stroke		
Stage up	Available*4			
Tube body	Binocular tube	Binocular S tube TC-T-TS (field number 22), Ergonomic ER tube TC-T-ER (field number 22)		
	Motorized eyepiece tube base unit for external PH (Ti2-T-BP-E)	Camera port (field number 16), Motorized PH turret with 4 motorized positions	—	
	Assist eyepiece tube base unit (Ti2-T-BA)	Assist camera (field number 22), Status detection		—
	Eyepiece tube base unit with port (Ti2-T-BC)	Camera port (field number 16)	—	Camera port (field number 16)
Transmitted illumination	Pillar for transmitted illumination (Ti2-D-PD)	Condenser vertical stroke: 66mm, Backward tilting up to 25 degrees, With field diagram and refocusing mechanism 2 filter slot positions (4 filter position option is also available with Filter Slider for transmitted illumination (Ti2-D-SF))		
	LED Lamphouse for dia illumination (Ti2-D-LHLED)	High power LED		
	Precentered Lamphouse (D-LH/LC)	100W halogen bulb (pre-centered)		
Condenser	Motorized condenser turret (Ti2-C-TC-E)	7 motorized positions (Ø37mm x4, Ø39mm x3), LWD/ELWD/CLWD/NAMC condenser lenses are supported	—	
	Intelligent condenser turret (Ti2-C-TC-I)	7 manual positions (Ø37mm x4, Ø39mm x3), Status detection, LWD/ELWD/CLWD/NAMC condenser lenses are supported	—	
	Condenser turret (TC-C-TC)	7 manual positions (Ø37mm x4, Ø39mm x3), LWD/ELWD/CLWD/NAMC condenser lenses are supported		
	ELWD-S condenser turret (TE-C)	4 manual positions, With ELWD condenser lens (NA0.3/OD65)		
	HNA condenser slider (Ti2-C-SCH)	2 manual positions (Ø37mm x1, Ø39mm x1), HNA dry lens/HNA oil lens are supported		
	Condenser lens	LWD (W.D.=30mm, NA=0.52), ELWD (W.D.=75mm, NA=0.3), CLWD (W.D.=13mm, NA=0.72), HNA dry (W.D.=5mm, NA=0.85), HNA oil (W.D.=1.9mm, NA=1.3), NAMC (W.D.=44mm, NA=0.4)		
Stage	Motorized stage (Ti2-S-SE-E, Ti2-S-SS-E)	Stroke X: ±57mm, Stroke Y: ±36.5mm, Max drive speed: approx. 25mm/sec, Magnetic sample holder	—	
	Stage (TC-S-SR, TC-S-SRF)	Stroke X: ±57mm, Stroke Y: ±36.5mm, Adjustable stroke range (3 levels) with adjusting pin, Long/middle/short handle options available		
	Gliding stage (TC-S-GS)	Stroke Ø20mm		
Nosepiece	Perfect Focus Unit with motorized nosepiece for Auto Correction Collar (Ti2-N-NDA-P)	5 motorized positions, Simple waterproof structure		—
	Motorized DIC sextuple nosepiece (Ti2-N-ND-E) Perfect Focus Unit with motorized nosepiece (Ti2-N-ND-P) Perfect Focus Unit with motorized nosepiece for MP (Ti2-N-NDM-P)	6 motorized positions, Simple waterproof structure		—
	Intelligent DIC sextuple nosepiece (Ti2-N-ND-I)	6 manual positions, Status detection, Simple waterproof structure		—
	Sextuple nosepiece (Ti2-N-N), DIC sextuple nosepiece (Ti2-N-ND)	6 manual positions, Simple waterproof structure		
Epi-filter turret	Motorized epi filter turret (Ti2-F-FLT-E, Ti2-F-FLTH-E)	6 motorized positions, Motorized shutter		—
	Intelligent epi filter turret (Ti2-F-FLT-I)	6 manual positions, Manual shutter, Status detection*5		
Filter wheel/shutter	Motorized BA filter wheel (Ti2-P-FWB-E)	7 motorized positions, High speed mode: 50ms, Low vibration mode: 100ms (movement time between adjacent positions)		—
	Motorized shutter (NI-SH-E)*6	12ms to open/close		
Epi-fluorescence attachment	EPI-FL module (Ti2-LA-FL), EPI-FL module for large FOV (Ti2-LA-FLL)	Supports fiber illuminator; includes 2-position filter slider and aperture diaphragm		
	Simple EPI-FL attachment (Ti2-F-FLS)	Supports both fiber illuminator and lamp house; includes 3-position filter slider		
	Field stop slider	Circular (Ti2-F-FSC), rectangular (Ti2-F-FSR), square (Ti2-F-FSS) aperture options		
Control unit	Controller, display device	Stage joystick (Ti2-S-JS), Tablet	Tablet	—
	Controller for Ti2-E (Ti2-CTRE)	USB/LAN interface, I/O function		—
Operating environmental		Temperature: 0°C+40°C, Humidity: 60% RH max. (at +40°C, no condensation), Indoor use only		

Motorized accessories have a status detection function

\*1 Motorized model with a bottom port

\*2 Limitations apply based on objective and filter cube choice, stage-up configuration,  
and illumination module, etc.

\*3 Tube base units with a port cannot be used with Ti2-A

\*4 Stage up kit is required. Please contact Nikon.



\*5 Status detection cannot be used when attached to the Ti2-U

\*6 NI-SH-CON Controller for Motorized Shutter is required for use with the Ti2-A/Ti2-U

# Dimensional Diagram

## Ti2-E



Double layer configuration with an Epi-FL module and a FRAP module

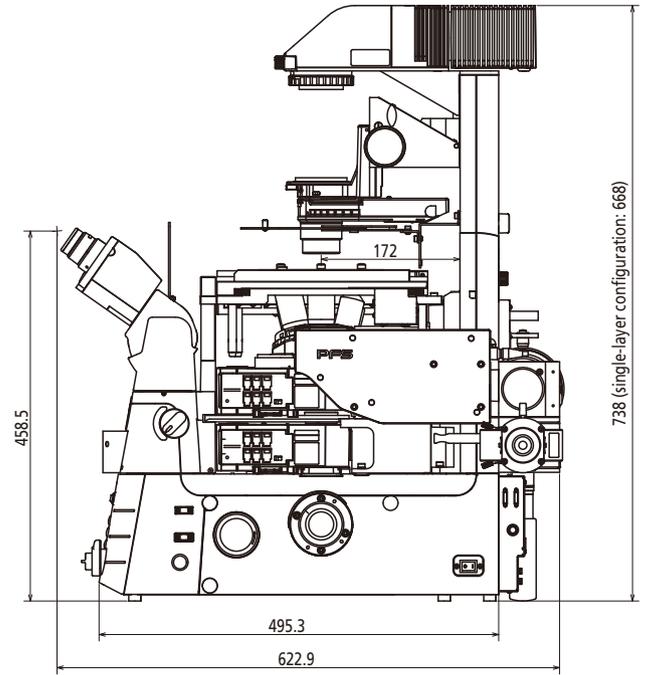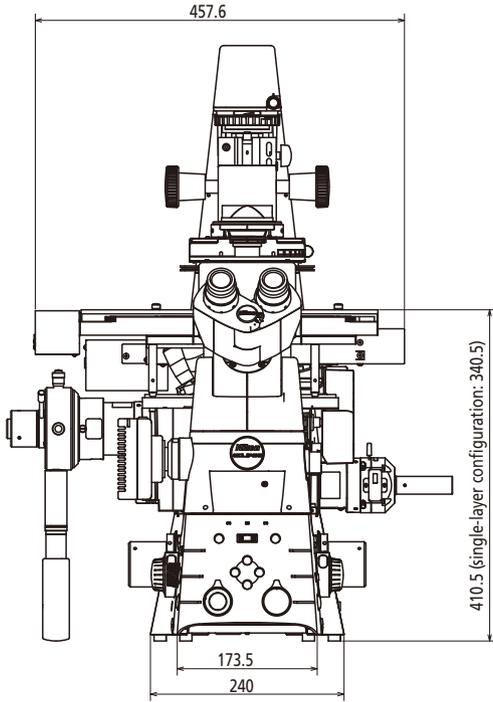

Unit: mm

## Ti2-A/U

Single layer configuration with an Epi-FL module

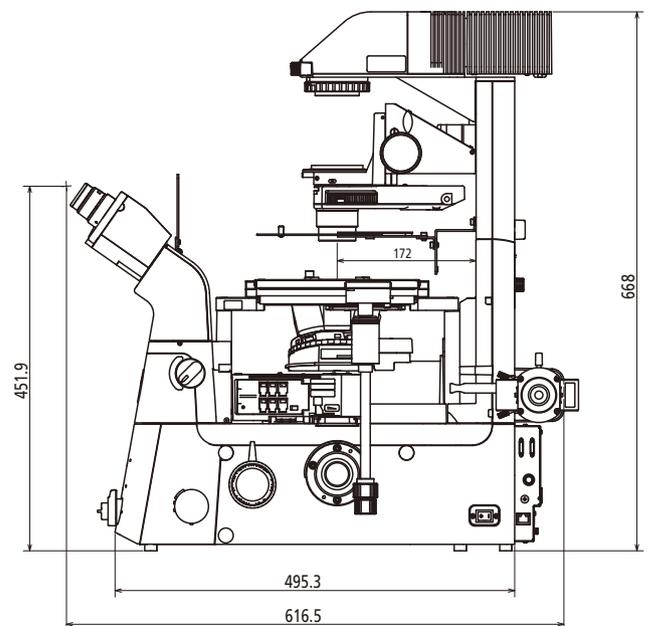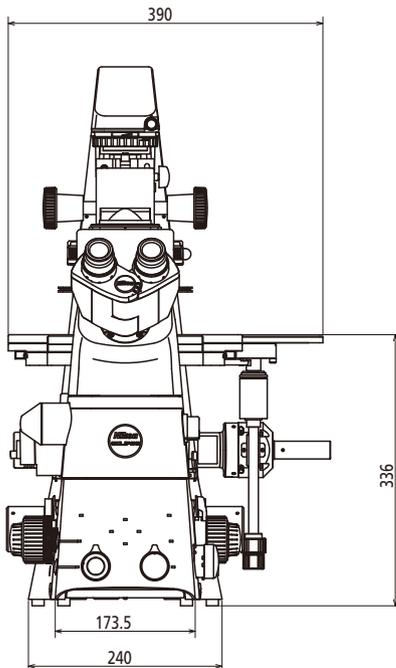

Unit: mm

Illustration is of Ti2-A

Enter the "Microscopy University" on the web  
and discover a whole new world. <http://www.microscopyu.com>

Specifications and equipment are subject to change without any notice or obligation  
on the part of the manufacturer. November 2016 © 2016 NIKON CORPORATION



**WARNING**

TO ENSURE CORRECT USAGE, READ THE CORRESPONDING  
MANUALS CAREFULLY BEFORE USING YOUR EQUIPMENT.

Monitor images are simulated.

Company names and product names appearing in this brochure are their registered trademarks or trademarks.  
N.B. Export of the products\* in this brochure is controlled under the Japanese Foreign Exchange and Foreign  
Trade Law. Appropriate export procedure shall be required in case of export from Japan.

\*Products: Hardware and its technical information (including software)



**NIKON CORPORATION**

Shinagawa Intercity Tower C, 2-15-3, Konan, Minato-ku, Tokyo 108-6290, Japan  
phone: +81-3-6433-3705 fax: +81-3-6433-3785  
<http://www.nikon.com/products/microscope-solutions/>

ISO 14001 Certified  
for NIKON CORPORATION

ISO 9001 Certified  
for NIKON CORPORATION  
Microscope Solutions Business Unit  
Industrial Metrology Business Unit

**NIKON INSTRUMENTS INC.**

1300 Walt Whitman Road, Melville, N.Y. 11747-3064, U.S.A.  
phone: +1-631-547-8500; +1-800-52-NIKON (within the U.S.A. only)  
fax: +1-631-547-0306

<http://www.nikoninstruments.com/>

**NIKON INSTRUMENTS EUROPE B.V.**

Tripolis 100, Burgerweeshuispad 101, 1076 ER Amsterdam, The Netherlands  
phone: +31-20-7099-000 fax: +31-20-7099-298

<http://www.nikoninstruments.eu/>

**NIKON INSTRUMENTS (SHANGHAI) CO., LTD.**

CHINA phone: +86-21-6841-2050 fax: +86-21-6841-2060  
(Beijing branch) phone: +86-10-5831-2028 fax: +86-10-5831-2026  
(Guangzhou branch) phone: +86-20-3882-0550 fax: +86-20-3882-0580

**NIKON SINGAPORE PTE LTD**

SINGAPORE phone: +65-6559-3651 fax: +65-6559-3668

**NIKON INSTRUMENTS KOREA CO., LTD.**

KOREA phone: +82-2-2186-8400 fax: +82-2-555-4415

**NIKON CANADA INC.**

CANADA phone: +1-905-602-9676 fax: +1-905-602-9953

**NIKON FRANCE S.A.S.**

FRANCE phone: +33-1-4516-45-16 fax: +33-1-4516-45-55

**NIKON GMBH**

GERMANY phone: +49-211-941-42-20 fax: +49-211-941-43-22

**NIKON INSTRUMENTS S.p.A.**

ITALY phone: +39-55-300-96-01 fax: +39-55-30-09-93

**NIKON AG**

SWITZERLAND phone: +41-43-277-28-67 fax: +41-43-277-28-61

**NIKON UK LTD.**

UNITED KINGDOM phone: +44-208-247-1717 fax: +44-208-541-4584

**NIKON GMBH AUSTRIA**

AUSTRIA phone: +43-1-972-6111-00 fax: +43-1-972-6111-40

**NIKON BELUX**

BELGIUM phone: +32-2-705-56-65 fax: +32-2-726-66-45



Nikon promotes the use of eco-glass  
that is free of toxic materials such as  
lead and arsenic.