

Place EasyClear where you want to install it.

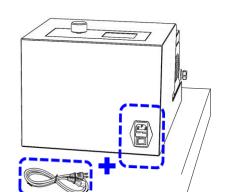


other end into a power outlet.


And Turn on the power switch.

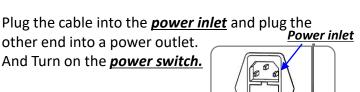

**Power Switch** 

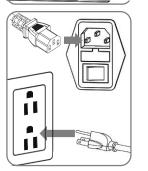

EasyClear Quick Guide

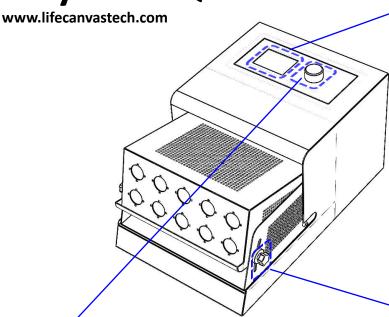


Press the **Temp.** button to set up the temperature and wait until the temperature stabilizes.

> 1 Press <u>Temp.</u> button 29.9 29.9 29.9 that matched with the tube hole. 2 Input target temperature to set-up 3 Enter to confirm


> > 30.0 29.9

- 4 Iterate 1~3 for each button for individual temp control
  - 5 Wait until A=B A .Current Temp.
  - B. Set-up Temp.


Regulate the shaking speed of the sample by adjusting the 'Shaking' knob and wait until the sample is processed.

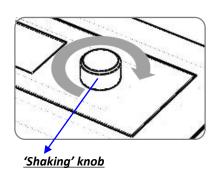

Place your sample into the tube and insert it into the device.

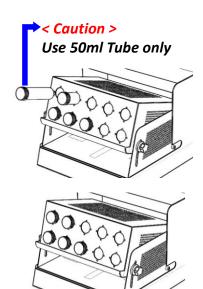

Regulate **the knob** on each side to adjust the height and the angle\*.

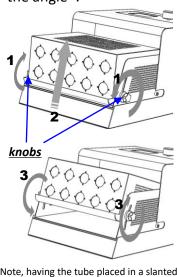

\*Note, having the tube placed in a slanted angle will help you use less reagents